

## OSHA Glove Selection Chart

OSHA Personal Protective Equipment Table 4 - <http://www.osha.gov/Publications/osh3151.pdf>

### Chemical and Liquid-Resistant Gloves

Chemical-resistant gloves are made with different kinds of rubber: natural, butyl, neoprene, nitrile and fluorocarbon (viton); or various kinds of plastic: polyvinyl chloride (PVC), polyvinyl alcohol and polyethylene. These materials can be blended or laminated for better performance. As a general rule, the thicker the glove material, the greater the chemical resistance but thick gloves may impair grip and dexterity, having a negative impact on safety. Some examples of chemical-resistant gloves include:

- **Butyl gloves** are made of a synthetic rubber and protect against a wide variety of chemicals, such as peroxide, rocket fuels, highly corrosive acids (nitric acid, sulfuric acid, hydrofluoric acid and red-fuming nitric acid), strong bases, alcohols, aldehydes, ketones, esters and nitrocompounds. Butyl gloves also resist oxidation, ozone corrosion and abrasion, and remain flexible at low temperatures. Butyl rubber does not perform well with aliphatic and aromatic hydrocarbons and halogenated solvents.
- **Natural (latex) rubber gloves** are comfortable to wear, which makes them a popular general-purpose glove. They feature outstanding tensile strength, elasticity and temperature resistance. In addition to resisting abrasions caused by grinding and polishing, these gloves protect workers' hands from most water solutions of acids, alkalis, salts and ketones. Latex gloves have caused allergic reactions in some individuals and may not be appropriate for all employees. Hypoallergenic gloves, glove liners and powderless gloves are possible alternatives for workers who are allergic to latex gloves.
- **Neoprene gloves** are made of synthetic rubber and offer good pliability, finger dexterity, high density and tear resistance. They protect against hydraulic fluids, gasoline, alcohols, organic acids and alkalis. They generally have chemical and wear resistance properties superior to those made of natural rubber.
- **Nitrile gloves** are made of a copolymer and provide protection from chlorinated solvents such as trichloroethylene and perchloroethylene. Although intended for jobs requiring dexterity and sensitivity, nitrile gloves stand up to heavy use even after prolonged exposure to substances that cause other gloves to deteriorate. They offer protection when working with oils, greases, acids, caustics and alcohols but are generally not recommended for use with strong oxidizing agents, aromatic solvents, ketones and acetates.

The following table from the U.S. Department of Energy (Occupational Safety and Health Technical Reference Manual) rates various gloves as being protective against specific chemicals and will help you select the most appropriate gloves to protect your employees. The ratings are abbreviated as follows: VG: Very Good; G: Good; F: Fair; P: Poor (not recommended). Chemicals marked with an asterisk (\*) are for limited service.

*Limited service	VG= Very Good	G= Good	F=Fair	P=Poor (not recommended)	
Chemical	Neoprene	Natural Latex or Rubber	Butyl	Nitrile Latex	
*Acetaldehyde	VG	G	VG	G	

Acetic acid	VG	VG	VG	VG
*Acetone	G	VG	VG	P
Ammonium hydroxide	VG	VG	VG	VG
*Amyl acetate	F	P	F	P
Aniline	G	F	F	P
*Benzaldehyde	F	F	G	G
*Benzene	F	F	F	P
Butyl acetate	G	F	F	P
Butyl alcohol	VG	VG	VG	VG
Carbon disulfide	F	F	F	F
*Carbon tetrachloride	F	P	P	G
Castor oil	F	P	F	VG
*Chlorobenzene	F	P	F	P
*Chloroform	G	P	P	P
Chloronaphthalene	F	P	F	F
Chromic Acid (50%)	F	P	F	F
Citric acid (10%)	VG	VG	VG	VG
Cyclohexanol	G	F	G	VG
*Dibutyl phthalate	G	P	G	G
Diesel fuel	G	P	P	VG
Diisobutyl ketone	P	F	G	P
Dimethylformamide	F	F	G	G
Diocetyl phthalate	G	P	F	VG
Dioxane	VG	G	G	G
Epoxy resins, dry	VG	VG	VG	VG
*Ethyl acetate	G	F	G	F
Ethyl alcohol	VG	VG	VG	VG
Ethyl ether	VG	G	VG	G

*Ethylene dichloride	F	P	F	P
Ethylene glycol	VG	VG	VG	VG
Formaldehyde	VG	VG	VG	VG
<b>Chemical</b>	<b>Neoprene</b>	<b>Natural Latex or Rubber</b>	<b>Butyl</b>	<b>Nitrile</b>
Formic acid	VG	VG	VG	VG
Freon 11	G	P	F	G
Freon 12	G	P	F	G
Freon 21	G	P	F	G
Freon 22	G	P	F	G
*Furfural	G	G	G	G
Gasoline, leaded	G	P	F	VG
Gasoline, unleaded	G	P	F	VG
Glycerine	VG	VG	VG	VG
Hexane	F	P	P	G
Hydrochloric acid	VG	G	G	G
Hydrofluoric acid (48%)	VG	G	G	G
Hydrogen peroxide (30%)	G	G	G	G
Hydroquinone	G	G	G	F
Isooctane	F	P	P	VG
Isopropyl alcohol	VG	VG	VG	VG
Kerosene	VG	F	F	VG
Ketones	G	VG	VG	P
Lacquer thinners	G	F	F	P
Lactic acid (85%)	VG	VG	VG	VG
Lauric acid (36%)	VG	F	VG	VG
Lineoleic acid	VG	P	F	G
Linseed oil	VG	P	F	VG
Maleic acid	VG	VG	VG	VG

Methyl alcohol	VG	VG	VG	VG
Methylamine	F	F	G	G
Methyl bromide	G	F	G	F
*Methyl chloride	P	P	P	P
*Methyl ethyl ketone	G	G	VG	P
*Methyl isobutyl ketone	F	F	VG	P
Methyl methacrylate	G	G	VG	F
Monoethanolamine	VG	G	VG	VG
Morpholine	VG	VG	VG	G
<b>Chemical</b>	<b>Neoprene</b>	<b>Natural Latex or Rubber</b>	<b>Butyl</b>	<b>Nitrile</b>
Naphthalene	G	F	F	G
Naphthas, aliphatic	VG	F	F	VG
Naphthas, aromatic	G	P	P	G
*Nitric acid	G	F	F	F
Nitromethane (95.5%)	F	P	F	F
Nitropropane (95.5%)	F	P	F	F
Octyl alcohol	VG	VG	VG	VG
Oleic acid	VG	F	G	VG
Oxalic acid	VG	VG	VG	VG
Palmitic acid	VG	VG	VG	VG
Perchloric acid (60%)	VG	F	G	G
Perchloroethylene	F	P	P	G
Petroleum distillates (naphtha)	G	P	P	VG
Phenol	VG	F	G	F
Phosphoric acid	VG	G	VG	VG
Potassium hydroxide	VG	VG	VG	VG
Propyl acetate	G	F	G	F
Propyl alcohol	VG	VG	VG	VG

Propyl alcohol (iso)	VG	VG	VG	VG
Sodium hydroxide	VG	VG	VG	VG
Styrene	P	P	P	F
Stryene (100%)	P	P	P	F
Sulfuric acid	G	G	G	G
Tannic acid (65%)	VG	VG	VG	VG
Tetrahydrofuran	P	F	F	F
*Toluene	F	P	P	F
Toluene diisocyanate	F	G	G	F
*Trichloroethylene	F	F	P	G
Triethanolamine	VG	G	G	VG
Tung oil	VG	P	F	VG
Turpentine	G	F	F	VG
*Xylene	P	P	P	F
<b>*Limited service</b>	<b>VG= Very Good</b>	<b>G= Good</b>	<b>F=Fair</b>	<b>P=Poor (not recommended)</b>